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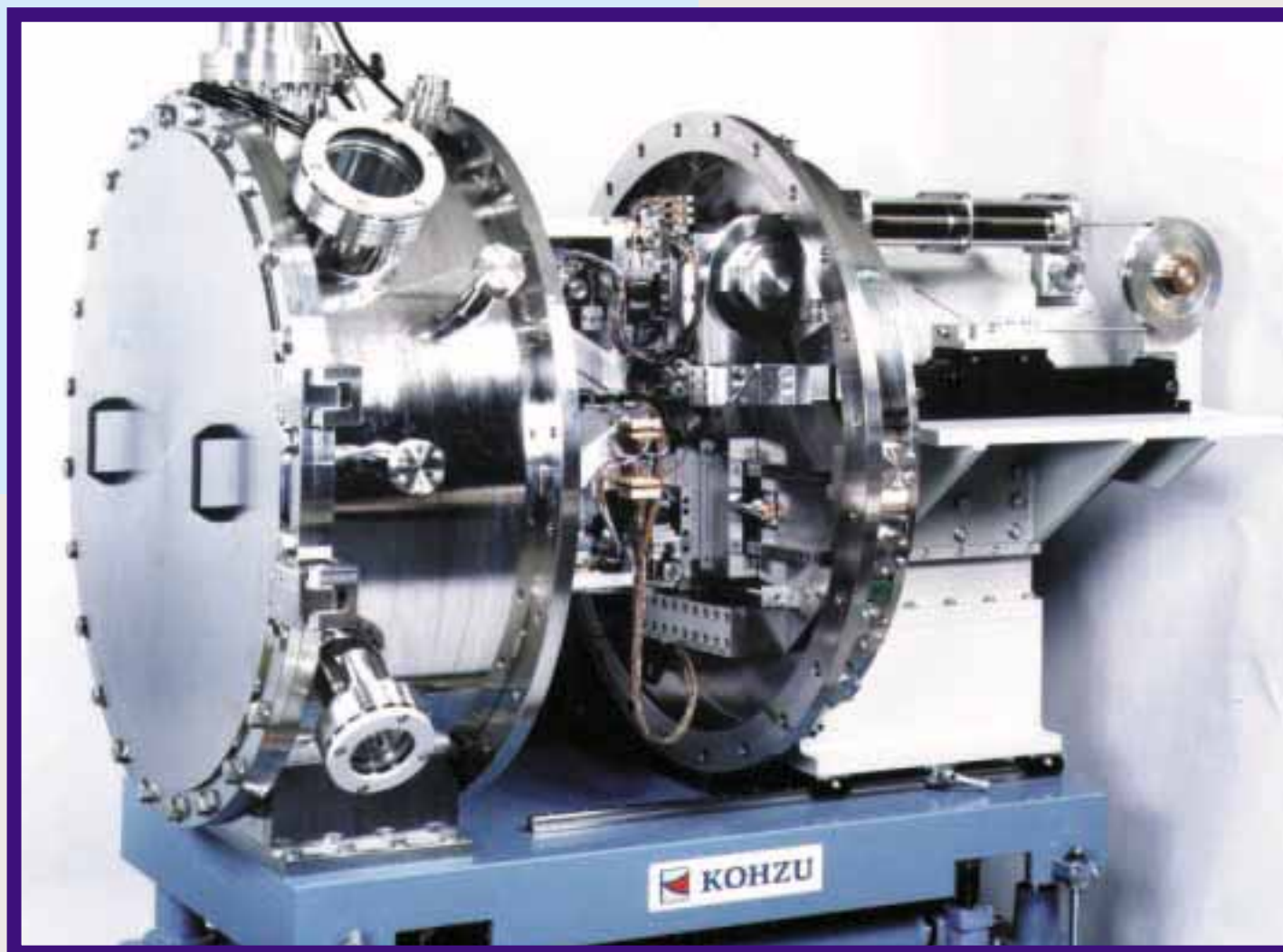
大型放射光施設 標準型2結晶X線分光器 STANDARD DOUBLE CRYSTAL X-RAY MONOCHROMATOR for SPring-8

General

As it is usual for synchrotron radiation facilities, X-ray double crystal monochromator produces a monochromatic exit beam, running parallel to incident white X-ray beam with fixed offset. It is required, that optical axis of the exit beam does not shift during a change of the exit beam energy produced by change of the crystal orientation, or the angle of incidence. In addition, it is desirable to change the energy of the exit beam by actuating least number of motors.

Here we offer our standard cam type monochromator with 1st and 2nd crystal holders placed on common rotating table. The design meets above described requirements and allows to have the exit beam in desired position. The monochromator is being produced in two modifications, either with "rotating inclined" or "variable inclined" geometry. Nevertheless, the design is the same except the crystal holder adjustment mechanics.

Main rotational axis coincides with the surface plane of the 2nd crystal. Shift of the 1st crystal position with respect to the incident beam, which would be caused by simple rotation of the table around the main axis is compensated by vertical motion of the 1st crystal guided by the cam. Then, the energy of the exit beam is determined only by the rotation angle without any displacement of the crystals with respect to the beam. Any thermally or mechanically caused distortions of the crystal parallelism can be corrected by precision tilting mechanism installed on the 1st crystal holder.

General Specification

Bragg angle range	3°~25°
Beam offset	30mm upwards
Vacuum degree	10 ⁻⁶ Torr

Every motor driven movement is secured by limiting switches.

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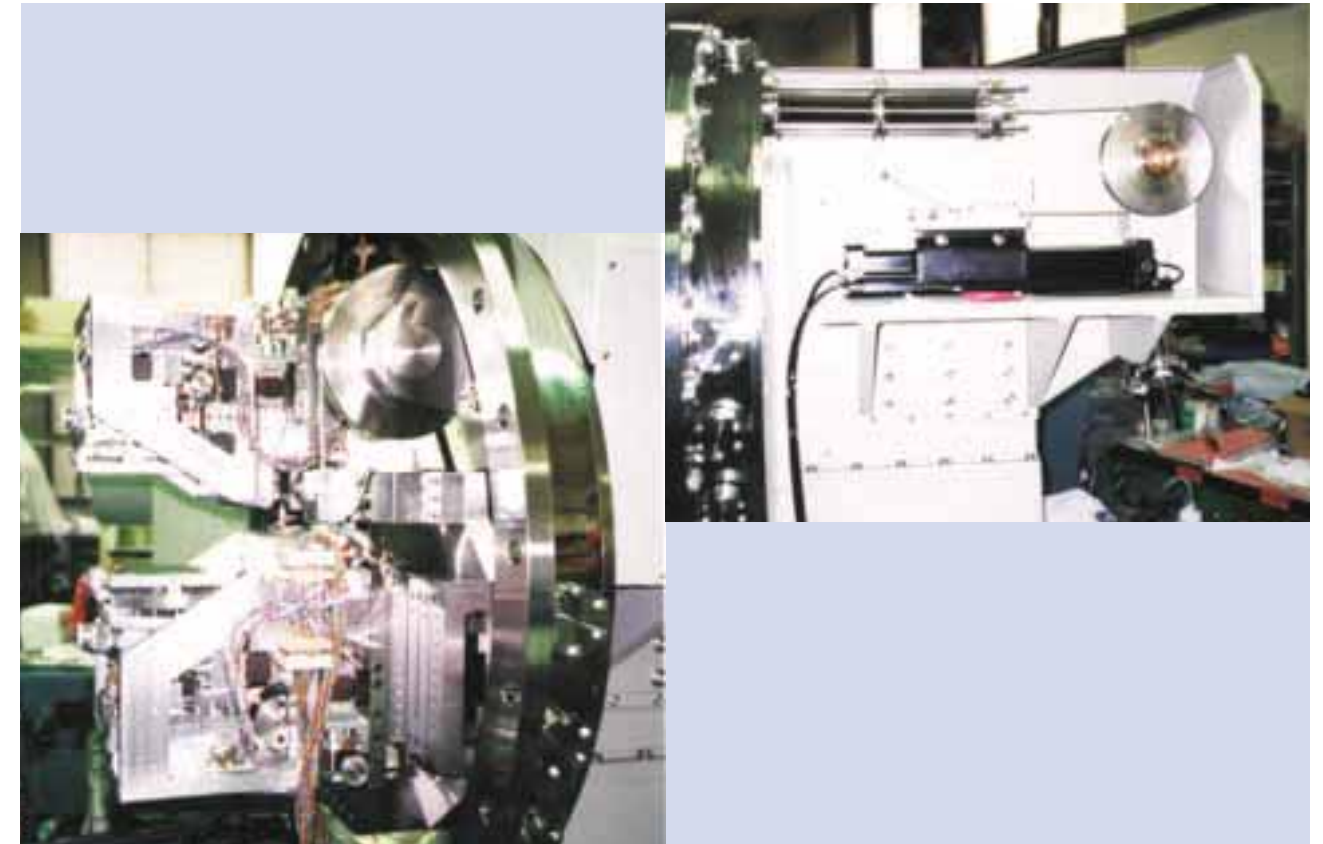
Main rotation axis is the most important part of the monochromator. It is because the high precision movement of the rotating table with all adjusting mechanisms in order to set the Bragg angle must be provided without breaking the vacuum. Nevertheless, our company possess the necessary technology, since we produced more than 30 monochromators based on the same principle, which are being successfully used* in Japan and abroad. Although the design of the monochromator is basically similar to that of other similar devices, it supposes the operation mainly at small angles of incidence. In this case, the distance between the crystals is large, and hence the load moment applied to the monochromator mechanism is large as well. In order to eliminate the resulting effect on the performance, the device is equipped with our peculiar moment load cancellation mechanism **.

* T.Ishikawa, SPring-8 Annual Report (JASRI, 1996) p.30.

** 特許申請中 patent application submitted


機構部仕様 Specification of the mechanism parts

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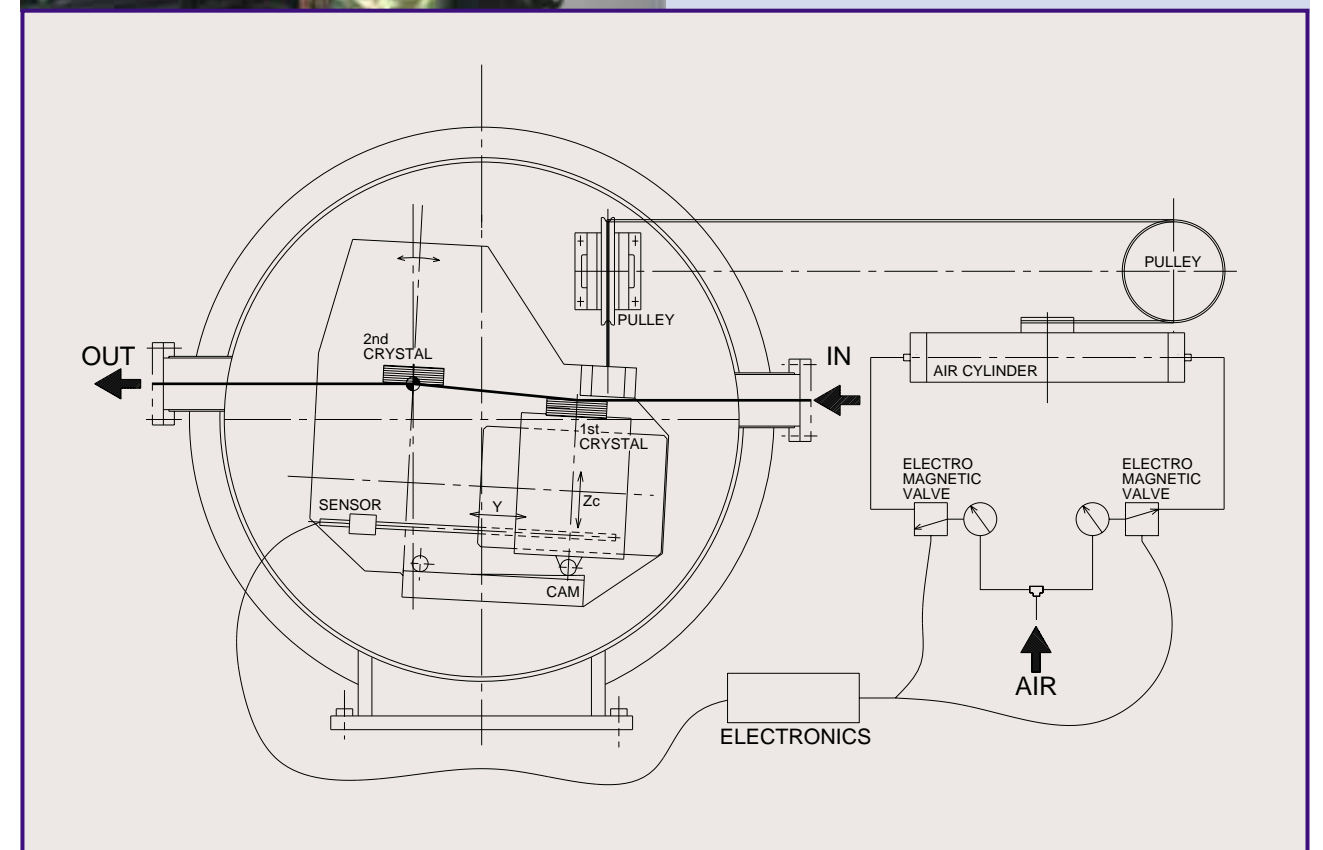- 1.In order to pass the cooling water tubes through the central axis, the doughnut-type encoder Heidenhain ERO-725 (1 reading) is being used.
- 2.The moment load cancellation mechanism is included.

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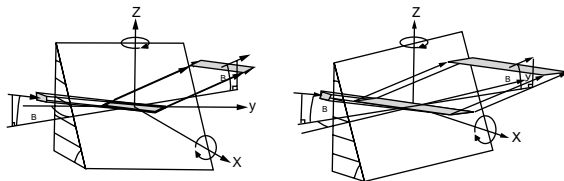
First crystal adjustment mechanism consists of the crystal holder adjustment mechanics, translation stage and the cam stage. The latter two parts secure desired position of the exit beam.


The holder adjustment mechanism, mounted on the horizontal 1-axis rotation stage, differs for the rotating inclined and variable inclined geometry except for x1-axis translation stage.

The translation stage motion can be controlled by vacuum compatible absolute linear motion magnetic sensor.

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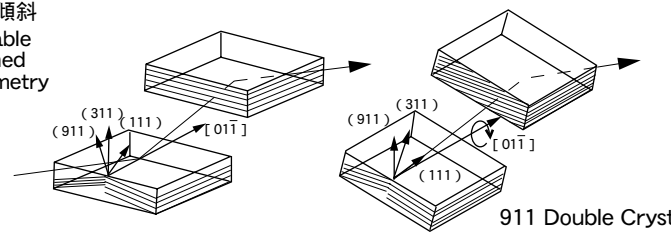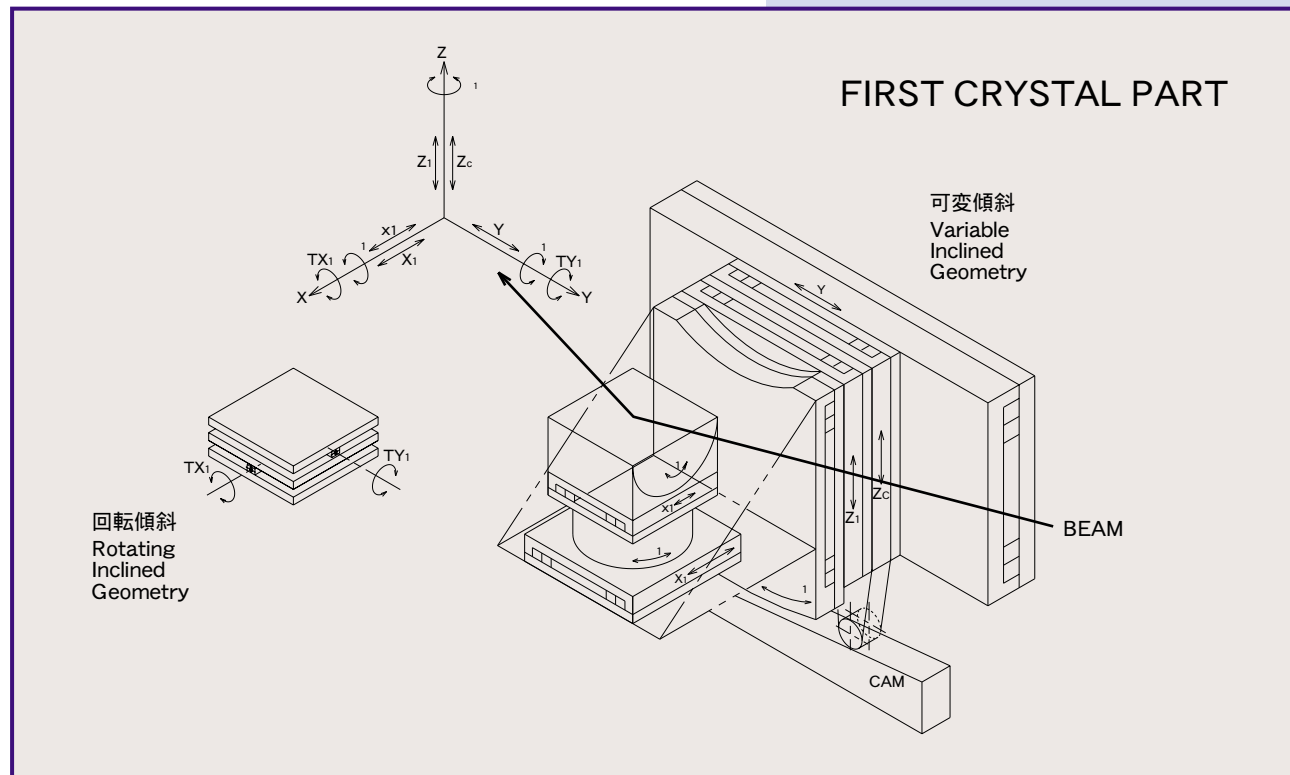
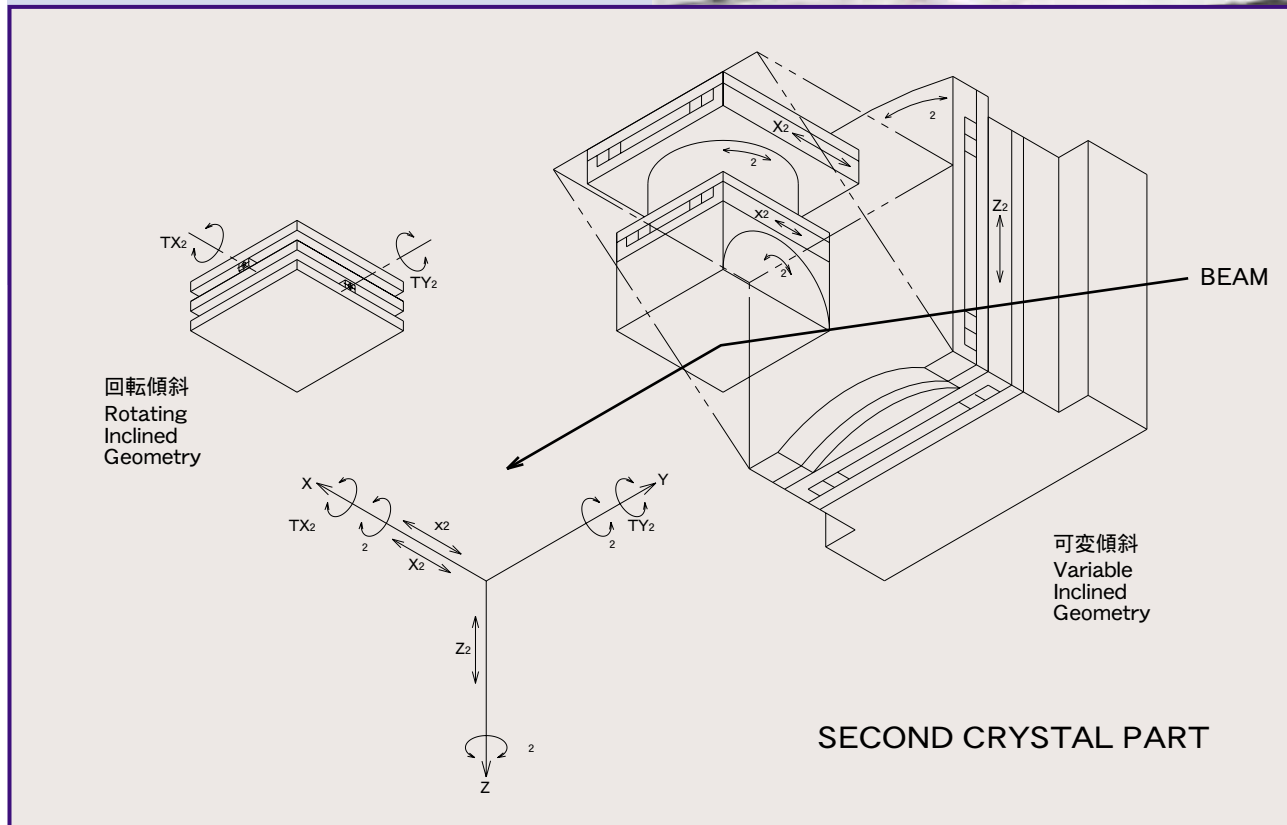
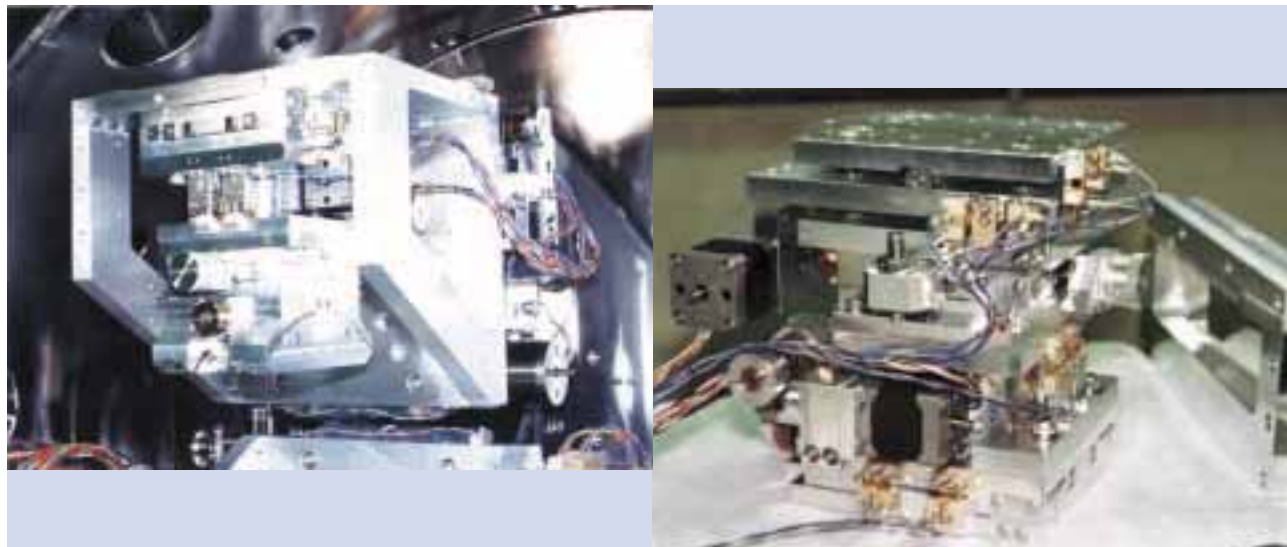
2nd crystal adjustment mechanism consists of the crystal holder motion stages. The function of this part is to adjust the relative position of the crystal surface with respect to the main rotation axis.

The holder adjustment mechanism, mounted on the horizontal 2-axis rotation stage, differs for the rotating inclined and variable inclined geometry except for x2-axis translation stage.

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4. 架台部 Supporting Structure

架台部構成 Supporting structure design

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Supporting structure is made of two parts. The upper part is a mild steel plate, which carries directly the vacuum chamber. The plate is equipped with rails to move the main rotation axis support with the mechanics.

The lower part is a welded structure support equipped with the Up/Down - Left/Right adjustment mechanisms.

調整軸仕様 Adjustment mechanics specification

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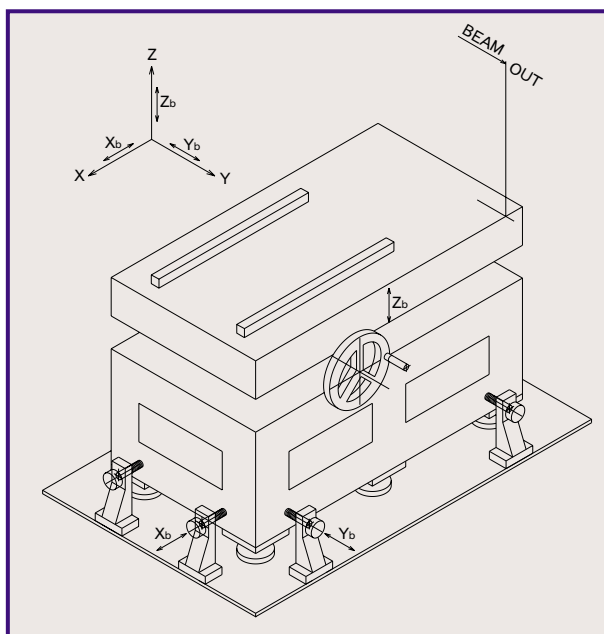
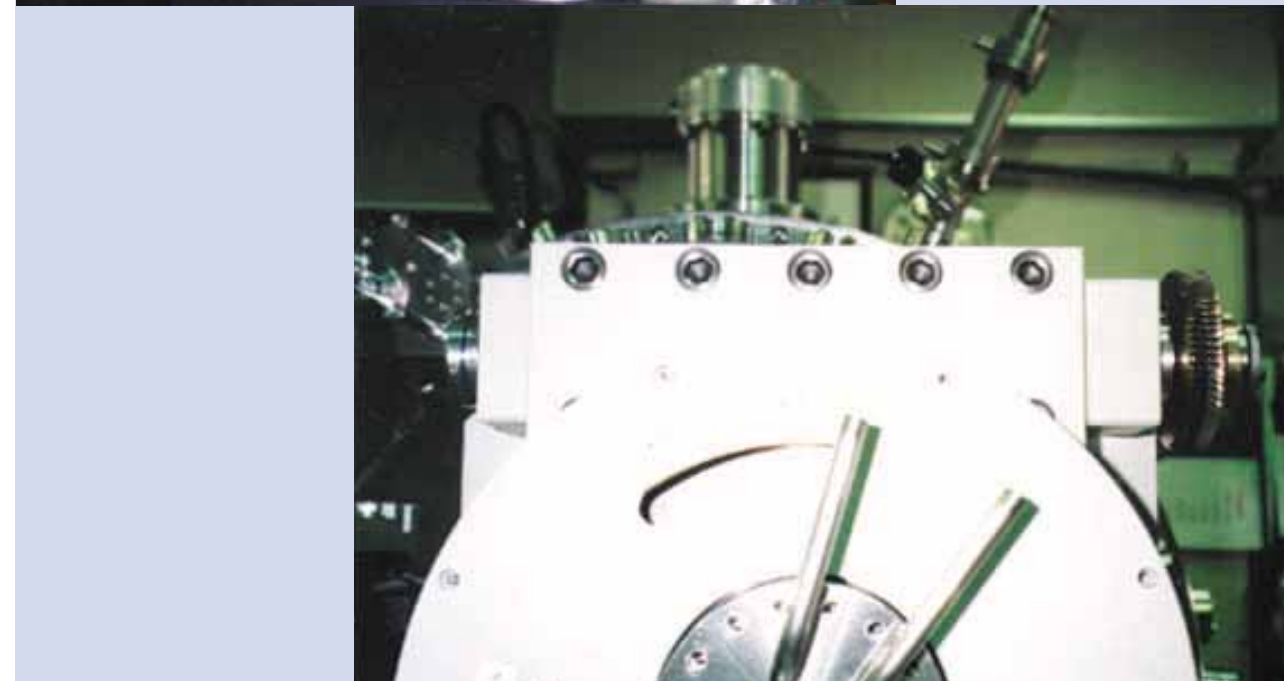
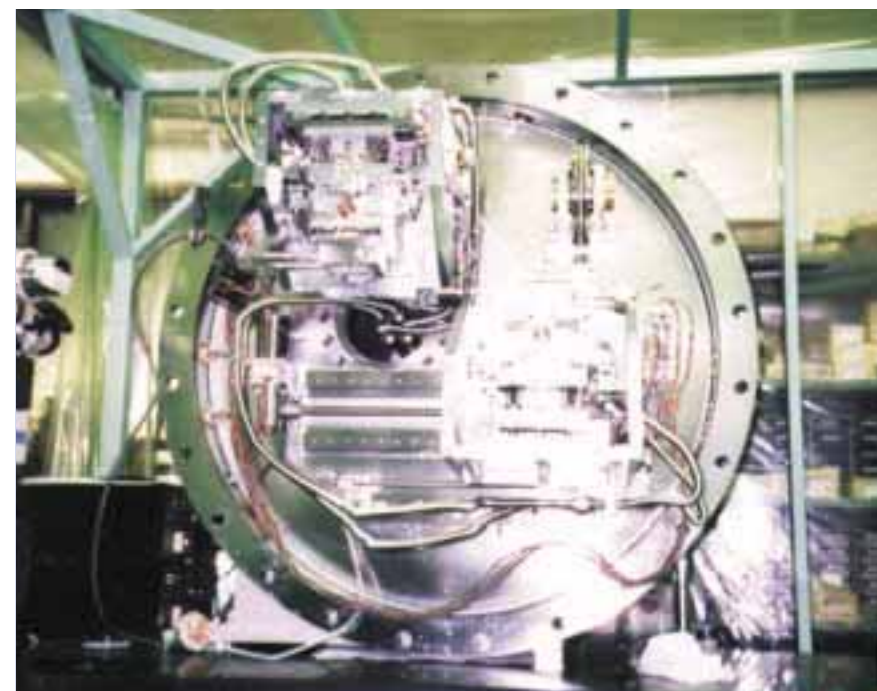
Flexible tubes are used inside the vacuum chamber for cooling water supply to the both crystals. The tubing connection in the vacuum is provided by VCR fitting. The joint size is 1/2 inch for the 1st and 1/4 inch for the 2nd crystal cooling system.

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Both crystal adjustment mechanisms are enclosed inside vacuum chamber body, covered with two large size flanges. Material of the chamber is SUS 304 stainless steel. The main axis shaft passes through the right side flange, if to look on the system from the beam source. The rotating table containing the adjustment mechanics of the both crystals is mounted on the vacuum side of the shaft.

The main axis shaft is passed to the vacuum chamber through a ferrofluid seal. Both left and right large size flanges are sealed by Viton seals.

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The monochromator is not baking type. Minimum achievable pressure is 10^{-6} Torr by turbomolecular pump.

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